

<b>Description</b>	<b>Precipitating polyclonal goat antiserum to rabbit immunoglobulins, IgG, IgA and IgM, heavy and light chains</b>	
<b>Product code</b>	GAR/Ig	
<b>Biological origin</b>	Goat	
<b>Physical form</b>	Delipidated, heat inactivated, lyophilized stable whole serum	
<b>Preservative</b>	No preservative added, as it may interfere with the antibody activity. No foreign protein added.	
<b>Immunogen</b>	Purified normal IgG and fractions containing IgA, and IgM isolated from rabbit serum. Freund's complete adjuvant is used in the first step of the immunization procedure.	
<b>Adsorption</b>	Immunoaffinity adsorbed using insolubilized antigens as required to eliminate antibody activity with other serum protein. The use of insolubilized adsorption antigens prevents the presence of excess adsorbent protein or immune complexes in the antiserum.	
<b>Identity &amp; Specificity</b>	The reactivity of the antiserum is directed to the major isotypes of the rabbit immunoglobulin system (classes and both light chain types) including antibodies to common determinants, to class and to the surface determinants of the common Fab portion, as tested by immunoelectrophoresis and double radial immunodiffusion against pooled serum and purified immunoglobulins.	
<b>Cross-reactivity</b>	Inter-species cross-reactivity is a normal feature of antibodies to immunoglobulins, since Ig of different species frequently share antigenic determinants. Cross-reactivity of this antiserum has not been tested in detail.	
<b>Protein concentration</b>	Total protein and IgG concentration in the antiserum are comparable to those of pooled goat serum. No foreign proteins added.	
<b>Antibody titre</b>	Precipitin titre not less than 1:32 when tested against normal rabbit serum in agar block titration.	
<b>Intended use</b>	In precipitating techniques as immunoelectrophoresis, single and double radial immunodiffusion (Mancini, Ouchterlony) to identify and measure immunoglobulins; as secondary antibody to precipitate the immunoglobulin in normal rabbit serum. This antiserum is not intended for use in non-precipitating antibody-binding or other highly sensitive assays. For such uses labelled and unlabelled cytochemical and immunoassay grade reagents meeting the necessary additional specificity and performance requirements are available.	
<b>Directions for use</b>	In immunoelectrophoresis use 2 µl or equivalent against 120 µl antiserum. In double radial immunodiffusion use a rosette arrangement with 10 µl antiserum in a 3 mm diameter centre well and 2 µl serum samples (neat and diluted) in 2 mm diameter peripheral wells.	
<b>Handling</b>	The lyophilized antiserum is shipped at ambient temperature and may be stored at +4°C; prolonged storage at or below -20°C. It is reconstituted by adding 1 ml sterile distilled water. Working dilutions are prepared by adding sterile phosphate buffered saline (PBS, pH 7.2). Repeated thawing and freezing should be avoided. Working dilutions should be stored at +4°C, not refrozen, and preferably used the same day. If a slight precipitation occurs upon storage, this should be removed by centrifugation. It will not affect the performance of the product.	
<b>Packing</b>	Vial with 1 ml lyophilized antiserum.	
<b>Storage / shelf life</b>	Lyophilized at +4°C	at least 10 years
	reconstituted at or below -20°C	3-5 years
	reconstituted at +4°C	7 days
<b>Caution</b>	This antiserum should be handled by qualified persons only and appropriate precautions should be taken in its handling and disposal, and of all associated materials. For <i>in vitro</i> laboratory research purposes only.	

**NORDIC IMMUNOLOGICAL LABORATORIES**

Langendijk 25, 5652 AX Eindhoven, The Netherlands

Tel. +31 630 070 625, Fax: +31 402 920 069

E-mail: [info@nordiclabs.nl](mailto:info@nordiclabs.nl)[www.nordiclabs.nl](http://www.nordiclabs.nl)