

Description	Precipitating polyclonal rabbit antiserum to human immunoglobulins IgG, IgA, IgM, IgD and IgE, heavy and light chains	
Product code	RAHu/Ig(GAMDE)	
Biological origin	Rabbit	
Physical form	Delipidated, heat inactivated, lyophilized stable whole serum	
Preservative	No preservative added, as it may interfere with the antibody activity. No foreign protein added.	
Immunogen	Purified normal Ig fractions isolated from pooled human serum. Freund's complete adjuvant is used in the first step of the immunization procedure.	
Adsorption	Immunoaffinity adsorbed using insolubilized Ig-depleted human serum fractions as required to eliminate antibody activity with other serum protein. The use of insolubilized adsorption antigens prevents the presence of excess adsorbent protein or immune complexes in the antiserum.	
Identity & Specificity	The reactivity of the antiserum is directed to the heavy and light chains of the immunoglobulin isotypes. It does not react with any non-Ig protein in human serum, as tested by immunoelectrophoresis and double radial immunodiffusion.	
Cross-reactivity	Inter-species cross-reactivity is a normal feature of antibodies to immunoglobulins, since Ig of different species frequently share antigenic determinants. Cross-reactivity of this antiserum has not been tested in detail.	
Protein concentration	Total protein and IgG concentration in the antiserum are comparable to those of pooled rabbit serum. No foreign proteins added.	
Antibody titre	Precipitin titre not less than 1:64 when tested against normal human serum in agar immunodiffusion block titration.	
Intended use	In immunoelectrophoresis and single and double radial immunodiffusion (Mancini, Ouchterlony) to identify the presence of immunoglobulins in human serum or other body fluids; to prepare an immunoabsorbent for the purification of human immunoglobulins from serum or plasma; as secondary antibody to precipitate the immunoglobulin in normal human serum. This antiserum is not intended for use in non-precipitating antibody-binding or other highly sensitive assays. For such use labelled and unlabelled cytochemical and immunoassay grade reagents meeting the necessary additional specificity and performance requirements are available.	
Directions for use	In immunoelectrophoresis use 2 µl or equivalent against 120 µl antiserum. In double radial immunodiffusion use a rosette arrangement with 10 µl antiserum in a 3 mm diameter centre well and 2 µl serum samples (neat and diluted) in 2 mm diameter peripheral wells.	
Handling	The lyophilized antiserum is shipped at ambient temperature and may be stored at +4°C; prolonged storage at or below -20°C. It is reconstituted by adding 1 ml sterile distilled water. Working dilutions are prepared by adding sterile phosphate buffered saline (PBS, pH 7.2). Repeated thawing and freezing should be avoided. Working dilutions should be stored at +4°C, not refrozen, and preferably used the same day. If a slight precipitation occurs upon storage, this should be removed by centrifugation. It will not affect the performance of the product.	
Packing	Vial with 1 ml lyophilized antiserum.	
Storage / shelf life	Lyophilized at +4°C reconstituted at or below -20°C reconstituted at +4°C	at least 10 years 3-5 years 7 days
Caution	This antiserum should be handled by qualified persons only and appropriate precautions should be taken in its handling and disposal, and of all associated materials. For <i>in vitro</i> laboratory research purposes only.	

NORDIC IMMUNOLOGICAL LABORATORIES

Langendijk 25, 5652 AX Eindhoven, The Netherlands

Tel. +31 630 070 625, Fax: +31 402 920 069

E-mail: info@nordiclabs.nlwww.nordiclabs.nl